



# OLYMPUS Stream 2.3.3

Release Notes  
September 2018

## Olympus Stream 2.3.3 Release Notes

ID	Title	Description
<b>R-2618</b>	Fix the problem that codebased activated licences expire after 30 days	Fix the problem with Stream 2.3/2.3.2 code based (by phone) activation procedure. Although there is no message during the code based activation process, all Material Solutions that are activated at the same time as the main license are actually NOT activated. This leads to an invalid Material Solution after 30 days of use. This problem does not affect Main licenses code based (by phone) activation but only Material Solutions. Alternative activation procedures are also perfectly working fine (online and by file exchange).
<b>R-2617</b>	Fix the problem that Windows 7-32 enters into an infinite boot loop	This release fixes the bug that appeared after April 4th 2018 when Microsoft has deployed a security update (KB4093118) for Windows 7-32 bit workstations in the world. We fix the issue that this security update create a conflict with the dongle installer (HASP driver) that is automatically installed with Stream (and cellSens) setup by removing the installation of the HASP driver on 32 bit operating systems.
<b>R-2622</b>	Fix the problem that report cannot be generated in Cast Iron Material Solution if several images from folder are used	The problem that a Cast Iron report containing multiple images from a folder generates only one image in the report is fixed.